



### Bottle/Vial/Syringe/Test Tube Washer

The Adelphi Bottle Washer is designed for rinsing the interior of containers with neck diameters between 9.5 mm and 100 mm, at a rate of up to 900 per hour. With the addition of optional nozzles, it can be used equally as effectively as a vial washer, syringe washer or test tube washer.

The unit consists of a spring loaded nozzle, which, when depressed by an inverted bottle, ejects jets of water which cleans every part of the inside of the container. The normal four armed rest is made from plastic and is easily cleaned or replaced.

Nozzles are screwed into the body and are interchangeable. The unit may be attached to the mains directly by plumbing (Model 3) or on a stand with a pipe connected to a tap (Model 2).

- A – part no. 6025000, Model 2 Bottle Washer
- B – part no. 6025100, Model 3 Bottle Washer, 0.5"-14 BSP threaded connection
- C – part no. 6025202, Sprinkler Nozzle
- D – part no. 6025201, Syringe Barrel & Test Tube Nozzle
- E – part no. 6025203, Vial Nozzle

